

TRINOCULAR STEREO MICROSCOPE

How to Assemble

The following are basic instructions on how to assemble the trinocular stereo microscope.




To Begin --


1.) Place the BASE on a strong, stable surface.




2.) Lower the BOOM ARM (with long side facing up) down onto the base POST.

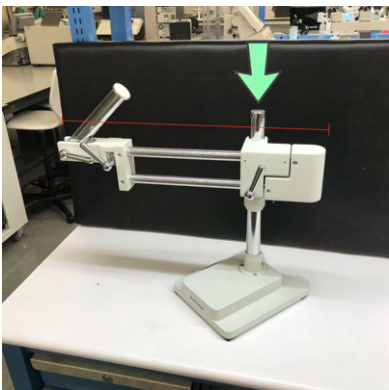


3.) Slide the BOOM ARM out to balance the assembly. Use the lever to secure it's position.

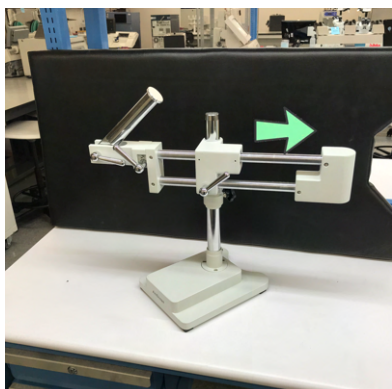


4.) Remove the threaded endcap from the PIVOT NECK and set aside temporarily.

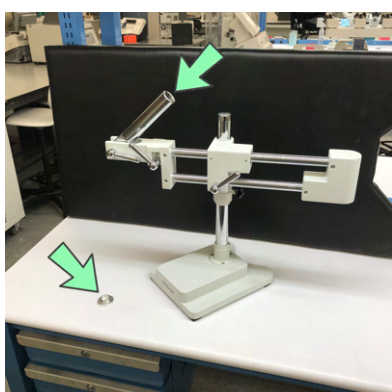

5.) Swing the PIVOT NECK into down position and tightly secure with lever.

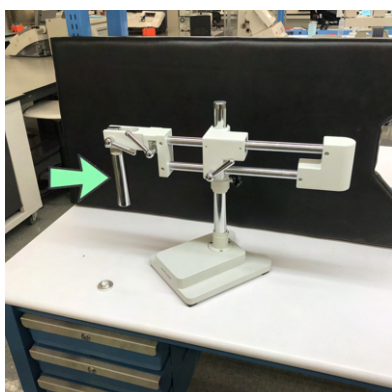

6.) Slide HEAD ASSEMBLY up onto NECK and screw endcap back into place at the end of the NECK.

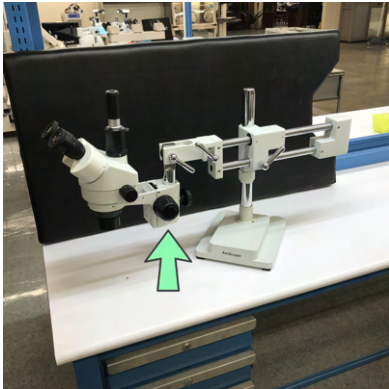

7.) Screw CAMERA onto center



8.) Slide the RING LIGHT up onto the lower part of the HEAD ASSEMBLY and secure it in place using the three thumbscrews.



[Click to HERE](#) to view item details on Alpha Props website